




LACTOSCAN MCCW

MILK ANALYZER Directly controlled by MS Windows Quick Start Guide

1 UNPACKING INSTRUCTIONS

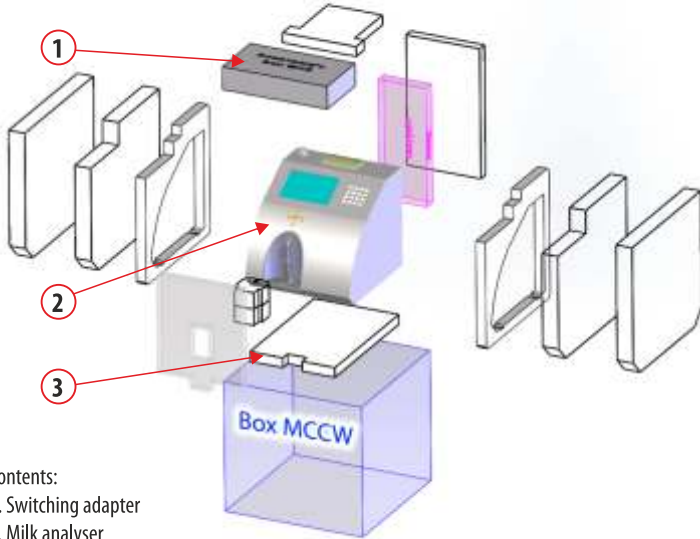

- Contents:
1. Switching adapter
 2. Milk analyser
Lactoscan MCCW
 3. Box with accessories

2 SAFETY INSTRUCTIONS

1. Read this manual carefully and make sure that you understand all the instructions.
2. For safety purposes the device is equipped with grounded power cable.
If there is no grounded electrical outlet where the device will be used, please, install such before using the device.
3. Place the device on leveled and stable plate.
In case it falls or is severely shocked it may be damaged.
4. Connect to the electrical network in such a way that the power cable to stay away from the side for accessing the device and not to be stepped on.
5. Do not disassemble the unit in order to avoid possible electrical shock.
In case of malfunction contact your local dealer.
6. Handle the liquids the device works with carefully, following all the instructions for their preparation.
7. Place the switching adaptor in such a way as to be protected from overflow and spillage of liquids.

3 ASSEMBLING THE DEVICE

Open the box. Carefully take out the contents of the box.
Place the machine on flat level surface

4 SWITCHING ON THE SYSTEM

1. Plug in the analyser power supply in one of the back panel inlets



Automatic cleaning (option)

2. Plug in any USB device which will be needed



3. Switch on the Lactoscan from the back



4. If you need to measure pH, follow the procedure below:

1. Loosen the nut anti-clockwise (1).
2. Pull up the stopper (2)
3. Carefully place the pH probe paying attention not to remove the sealing O-ring (3)
4. Place the probe with the nut in the hole (C & D) and tighten it.



If you'll not use the analyser more than 2 days, you must take out the probe and to place the stopper back.
The pH probe must be stored separately as per the instructions of point 2.3 Appendix 3.

5. Power On the Tablet
Wait until the software opens



6. If everything is connected properly you should receive a "Ready to Start" message on the screen



7. Fill the glass with a milk sample and place it in the analyser



8. Start a measurement by pressing the calibration name on the screen for the type of milk you are going to measure (if you are measuring cow milk, press the Cow Button)



9. When the measurement is started on the screen appears



Where: Position 1: The seconds left until the measurement is finished
 Position 2: Temperature of the sample
 Position 3: The calibration channel that is being used
 Position 4: Message about what calibration you are using for measurement

10. When the measurement is finished



Where: Position 1: Highlighted the last used calibration
 Position 2: Result collector for all the measurements done since the start of the program
 Position 3: Measure count since the start of the program
 Position 4: Archive file for the results
 Position 5: Fields for the parameter results of the last measurement

12. Continue work by placing a new sample in the glass and press the calibration buttons

13. When finished work put clean water in the glass and start cleaning by pressing the cleaning button



14. Please refer to the user manual about the proper way of cleaning the machine

IMPORTANT THE MAIN REASON FOR MALFUNCTIONING OF THE DEVICE IS THE BAD CLEANING OF THE SYSTEM AFTER MAKING ANALYSIS. In case of malfunction due to the bad cleaning of the analyser your guarantee is not valid anymore and any repair has to be paid.

5 TURNING OFF THE SYSTEM

1. To turn off the tablet select **File -> Shutdown**



2. To exit the program select **File->Exit** to close the software to use the tablet



3. Turn off the analyser from the back panel switch

*NOTE: If you have any questions about any of the procedures, please contact Milkotronic LTD

6 OPTIONS AND ACCESSORIES



7 MORE INFORMATION

You can download the Instruction Manual of the device from our website http://lactoscan.com/editor/uf/manuals/MCCW/Lactoscan_MCCW_Eng.pdf You can also find it in the Help Menu of the analyser:



The Instruction Manual of the Data Base Application can be downloaded from the Help Menu of the analyser.

8 SUPPORT

For any additional information and further question regarding the servicing of the device, contact us at: mainservice@lactoscan.com

HEAD OFFICE:
 4, Narodni Buditeli Str.
 8900 Nova Zagora
 BULGARIA
 Phone/Fax: + 359 457 67082

MILKOTRONIC LTD
office@lactoscan.com
www.lactoscan.com
www.milkotronic.com